



## Workspace Templates Pollen Viability

This template download collection contains measurement templates for pollen viability measurements of the following species:

- Artichoke
- Barley
- Broccoli
- Carrot
- Cauliflower
- Corn
- Cucumber
- Daisy
- Fennel
- Melon
- Onion
- Pepper
- Sorghum
- Squash
- Sunflower
- Tomato
- Watermelon
- Wheat
- Zucchini

For pollen counting, download the pollen counting collection from the website.

Our *Quick Guides* contain valuable tips for experimental setup.

Download Viability  
Template Collection

If a template for a specific species you might need is missing, please contact us at [support@amphasys.com](mailto:support@amphasys.com)

### Download at

[amphasys.com/download](https://amphasys.com/download)

### Download and use

- Download the template collection
- Unzip the template collection
- Open the template collection folder
- Create a copy of the desired template
- Rename the copy. This will be your workspace
- Open AmphaSoft and load the workspace

### Tips

- Use our Quick Guides *Getting Started with Pollen Analysis* and *Pollen Analyzer Startup Procedure* to get started
- Keep on copying the original template folder to create individual workspaces

### Support

If you need support for the template management and cannot find a solution in the *AmphaSoft User Guide* or our *Quick Guides*, contact us at

[support@amphasys.com](mailto:support@amphasys.com)