



New Large Pollen Kit For Ampha Z32 Pollen Analyzer

Measuring big pollen is now possible with the new 400 µm AmphaChip. Discover the new «Large Pollen Kit» for the Ampha Z32 Pollen Analyzer and measure big pollen with a diameter between 100 µm – 200 µm.

Product no / name	11.954 / Large Pollen Kit
Species recommended	<ul style="list-style-type: none"> ▪ Pumpkin, Squash, Zucchini (Cucurbita maxima, C. pepo, C. moschata and other C. species) ▪ Cotton, Ocra, Hibiscus and other species of Malvaceae family
Pollen size	From 100 µm up to 200 µm
Kit content	2 AmphaChips 400 µm 50 filters, 300 µm 50 filter adapters 2 bottles of AF 8 or AF 9 buffers
Buffer	Works only with new AF Buffer 8 and 9
Technical details of chip	Channel size 400 µm width x 300 µm height
AmphaSoft version required	AmphaSoft 2.1.6
Additional information	Species specific templates needed for measurement, templates can be downloaded from Amphasys website: www.amphasys.com/downloads
Can be ordered as of	Immediately, please contact: order@amphasys.com
Can be delivered as of	January 2020



Download and use

- [Download](#) the template collection
- Unzip the template collection
- Open the template collection folder
- Create a copy of the desired template
- Rename the copy. This will be your workspace
- Open AmphaSoft and load the workspace

Tips

- Use our [Quick Guides](#) *Getting Started with Pollen Analysis* and *Pollen Analyzer Startup Procedure* to get started
- Keep on copying the original template folder to create individual workspaces

Price	
List price as of 1.1.2020	CHF 1'465.00